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The battery box 251580 is used for powering Frederiksen's microphones and other analog sensors when they are connected to equipment that doesn't provide power. A typical example is connecting a microphone (248600 or 251570) to an oscilloscope or an AC voltmeter.

The sensor input is a 6-pin modular socket. We have included a flat cable with modular plugs which fits e.g. the microphones mentioned.

Older microphones and sensors with DIN plugs cannot be used with 251580 but connects to the older version of the battery box 251560.

Use

The battery box 251580 turns on automatically when connected through the modular socket. This is indicated by an LED.

Don't forget to remove the plug from the socket when the equipment is no longer used.

The analog signal is accessed through the two 4 mm safety sockets.

Battery change

The battery box uses a single 9 V battery (item No. 351010).

Remove the rubber pads at the bottom of the box. This exposes four small crosshead screws. Remove these and the bottom comes loose and the battery can be replaced.

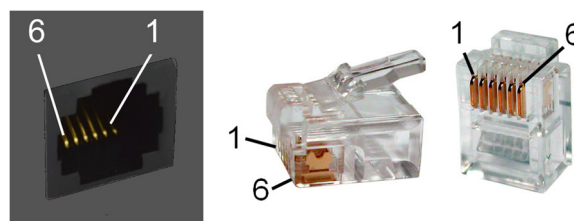
When reassembling the box, take care not to tighten the screws more than necessary. Nudge the rubber pads back in place again.

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© Frederiksen Scientific A/S